## **UM-CHOR-1 Growth Conditions**

## **Growth conditions:**

IMDM:RPMI 1640 mixed 4 parts IMDM:1 part RPMI 10% FBS 1X NEAA Pen-Strep (optional)

Cells are grown on regular plastic and are split 1:2, when nearing confluence, about ever 3-4 days. It is best to wait an extra day if you are not sure rather than split too early. Cell should never be split back further than 1:4.

## **Freezing Media:**

70% complete growth medium 20% fetal bovine serum 10% DMSO